

EUTHANEX INSTRUCTION (Blue Line)

1. Ensure green flowmeter for individual cages is turned off.
2. Remove cage lids (leave wiretop in place).
3. Place cages in Euthanex chamber & close.
4. Select Adults (15 minute cycle) or Neonates (60 minute cycle).
5. Press RUN.
6. Do not open door until cycle is finished.
7. Press STOP to abort cycle if necessary.

SC #237
Eff 2/19

3.5" x 5" horizontal template
Inner box dimensions 3.2" x 4.6"
Outer box dimensions 3.5" x 4.8"

EUTHANEX INSTRUCTION (Green Line)

1. Ensure green flowmeter for individual cages is turned off.
2. Remove filters from the Green Line lids
3. Place cages in the Euthanex chamber & close.
4. Select Adults (15 minute cycle) or Neonates (60 minute cycle).
5. Press RUN.
6. Do not open door until cycle is finished.
7. Press STOP to abort cycle if necessary.

SC #237
Eff 2/19