

SAFETY FEATURES

- **To prevent entrapment the buddy-system must be used.**
- **Two individuals must be present each time the bulk sterilizer is started.**
- **The inside of the autoclave is observed to confirm that no one is inside prior to activating.**
- **Both individuals will remain outside the autoclave to ensure that no one is inside as the door is closed and the sterilizer cycle is being activated.**
- **By design the door to the sterilizer only closes very slowly to reduce the possibility of entrapment.**
- **Emergency Shut-Off (red button) located on both sides of the autoclave.**
- **Emergency Electrical Cut-Off located to left of unit on prep-side.**

SC #259
Eff 7/19

SRB Large Bulk Sterilizer